

Important notes

The return transport - one call is enough!

The return transport is carried out by our logistics partner.

Please observe the export regulations of your country.

GO! Express & Logistics GmbH

Kunden-Nr: 103977

Phone Germany: 0800 1112828

Phone abroad: +49 7131 155650

Please note on the day of sampling:

- Registration of the collection until 12 o'clock at the latest
- Collection takes place until 2 pm
- Last possible return day is always Thursday
- No collection before holidays

Content:

- 20 Rodac plates for the examination of surfaces
- 1 waterproof pen
- 1 Bottle of hand sanitizer
- 1 laundry net with temperature logger
- 1 Cooling element

Do you have any questions:

Write to us at:
info@hymo-box.de



Weitere Informationen
finden Sie unter:
www.hymo-box.de

Temperature check Washing process

Hygiene monitoring made easy

The
**HyMo-
Box**

Take your hygiene management into your own hands with the HyMo-Box!

Using temperature data loggers, you can easily check your washing processes in practical operation and document the proper condition of your machines.

With the additional swab tests, you can verify the machine, the wash ware or the environment for the low germ count after the washing process. Our accredited laboratory provides you with the germ counts of your individually defined sampling points.

Please note the instructions for return transport on the reverse side.



Step by step guide

for the HyMo-Box temperature check

Your advantages: You can easily and reliably monitor your cleaning and disinfection measures. It's all inclusive - there are no hidden costs for you. Our DAKkS accreditation guarantees you the highest quality of our services.



- Open HyMo-Box.
- Place cooling element in the refrigerator.
- In case of later processing also store the box with the RODAC plates in the refrigerator.




- Define day of performance of tests and organize return transport for the same day.
- Attention: registration of the return transport must be made before 12 noon, the collection of the shipment will be made by 2 pm, no return transport on Fridays.



3

- Remove the laundry net from the packaging.
- Place in the standard washer-extractor or cycle washing machine. Start the washing program.

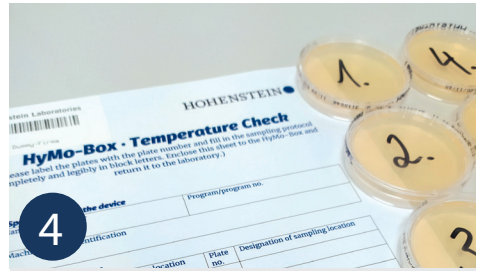

4

- Define sampling sites for the determination of the surface bacterial count and enter them in the sampling plan.
- Label RODAC plates with numbers according to the sampling plan.



5

- Disinfect hands.
- Open the lid of a RODAC plate.
- Press the culture medium surface for approx. 5 seconds under slight pressure on the area to be examined - do not move the RODAC plate if possible.
- Close the lid again.
- Proceed with all further RODAC plates according to your sampling plan.



6

- At the end of the washing program, remove the net from the laundry and place it in the plastic bag provided.
-



7

- Stack the RODAC plates back into the box. Place this in the box, place the laundry net in the cooler, close the box and place in the outer box.
- Place sampling plan on box and tape box closed.